


Vacuum oven VO with TwinDISPLAY AtmoCONTROL software

Model sizes:
29 / 49 / 101
+20 °C to +200 °C
5 mbar to 1100 mbar
Accessories: lower pump chamber and energy-efficient vacuum pump

**VACUUM OVEN VO** The high-performance turbo dryer impresses with its many intelligent Memmert features for gentle drying and precise, rapid temperature control: digital pressure control, directly heated and individually controllable thermoshelves, and simple programming via ControlCOCKPIT or AtmoCONTROL software. Combined together, the speed-controlled vacuum pump and the vacuum oven VO are an unbeatable energy-efficient pairing. The pump fits neatly inside the matching lower chamber.

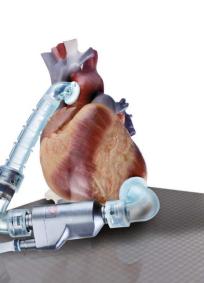






### Unique precision: Memmert VO direct heating

Available only from Memmert: multi-level sensing and heating. For really short heating-up and processing times, heating is provided via individually positionable thermoshelves with integrated shelf heating and sensors. The separate control circuits react precisely to different loads or humidity levels and ensure the setpoint temperature is consistently maintained. Due to the direct contact between the heating and the chamber load, there is practically no loss of heat. Each thermoshelf can be calibrated individually.

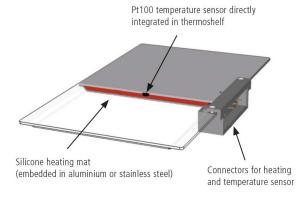

Removable thermoshelf with direct heating system and sensor

Multi-level sensing and heating

## Optional vacuum pump saves around 70 % energy


The speed-controlled chemically resistant Memmert vacuum pump is automatically detected by each vacuum oven VO. Thanks to intelligent speed control, it controls the setpoint with great precision. The energy efficiency is also obvious, with measurements showing energy savings of around 70 % in ramp mode compared with vacuum pumps that are not controlled; it is even possible to achieve higher savings at constant vacuum levels. The final vacuum level of up to 2 mbar favours a wide range of applications, while pump control (based on individual requirements) significantly extends the service life of membranes. If another vacuum pump or a central vacuum supply is connected, vacuum control is achieved via solenoid valves.



# Turbo drying thanks to vacuum cycles

Digitally controlled vacuum cycles, during which the working chamber is intermittently vented at short intervals, can achieve further significant reductions in drying times. The AtmoCONTROL software makes it quick and easy to program ramps with different temperature and vacuum setpoints.



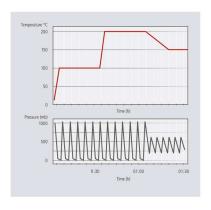

Example of ramp programming

## Convenience in a package: the Premium Module

The basic version of the vacuum oven VO features a thermoshelf and two thermoshelf connectors (VO29: 1 thermoshelf connector). The Premium Module includes the option for switching to inert gas, a programmable, digitally controlled gas inlet with flow reduction; there is also the MobileALERT option with separate error messages for temperature and pressure as well as (depending on the appliance size) additional thermoshelves and thermoshelf connectors (see the technical data for details).

#### **VACUUM OVENS VO**

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1)

Standard units are safety-approved and bear the test marks: C €







Interior:

Stainless steel interior, material 1.4404 (ASTM 316 L), hermetically welded, with removable mountings at the sides for cleaning, including thermoshelf guide bars, as well as mounting on top to avoid turbulences

Housing:


Textured stainless steel, rear zinc-plated steel, intuitively operated TwinDISPLAY (TFT colour displays) with touchscreen, safety glass door with inner bullet-proof glass and external anti-splinter screen

Mains cable with plug (German type) Connection:

Installation: 4 feet

Interfaces:



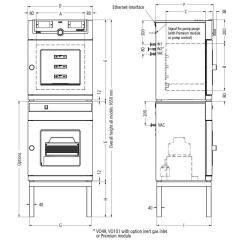

			or i remain module				
Model sizes/Descrip	ption			29	49	101	
Stainless steel interior	Volume	â	approx. I	29	49	101	
	Width	(A) mm		385		545	
	Height	(B)	mm	305	385	465	
	Depth	(C)	mm	250	330	400	
	Distance between thermoshelves		mm	75		95	
	Maximum load per oven	aj	pprox. kg	kg 40 6		0	
	Max. number of thermoshelves		number	1		2	
	Max. number of thermoshelves (with premium module)		number	2 4		4	
	Max. loading per thermoshelf		kg	20			
Textured stainless steel exterior	Width	(D)	mm	55	0	710	
	Height	(E)	mm	607	687	767	
	Depth (without door handle, depth of handle +38 mm)	(F)	mm	400	480	550	
	Safety glass door: Textured stainless steel frame with spring-loaded safety glass on inside and anti-splinter screen ESG on outside of door			•			
	Door Seal: Endless silicone profile seal				•		
Standard equipment	Thermoshelves – aluminium eloxadised , mat. 3.3547 (ASTM B209) – with integrated large-area heating including local temperature sensing (Pt100, 4-wire-circuit); individual overtemp. protection for each shelf. Further data see stainless steel number inner working chamber	İ	number	1			
	Works calibration certificate (measuring point in the middle of the individual shelf for +160 °C at 20 mbar pressure): a separate certificate is prepared for each thermoshelf ordered and shipped together with the vacuum oven		°C	•			
Temperature	Temperature sensors Pt100 Class A in 4-wire circuit individually for each thermoshelf				•		
	Working temperature range		°C	at least 5 above ambient temperature to +200			
	Setting temperature range		°C	+20 to +200			
	Setting accuracy		°C	up to 99.9: 0.1 / from 100: 0.			
	Temperature variation in time (aluminium thermoshelf)		K	$\leq \pm 0.3$			
	Temperature uniformity (surface) at +160 °C/20 mbar (aluminium thermoshelf)		K	≤ ± 2.5			
Pressure (vacuum)	Vacuum connection with small flange DN16, and gas inlet with fresh air supply				•		
	Digital electronic pressure control for a speed-controlled vacuum pump. Tubing for vacuum, air and inert gas are made of material 1.4571 (ASTM 316 Ti). Adjustable from 5 mbar up to 1100 mbar. Programmable, digitally controlled inlet for air				•		
	Pump control: optimised rinsing procedures for the pump membranes as well as signal output for pump ON/OFF				•		
	Rapid air intake for door opening without alteration of selected vacuum setpoint				•		
	Permitted final vacuum		mbar	0.01			
	Maximum leakage rate		bar/h	0.01			
Control technology	Digital over- and undertemperature monitor				•		
	Temperature monitoring band automatically linked to the setpoint (ASF)				•		
	Monitor relay for reliable heating cut-off in case of fault				•		
	Mechanical temperature limiter (TB)				•		
	Multi-Level-Overtemperature-Protection (MLOP) for each thermoshelf						

Model sizes/Des	•			29	49	101	
Further data	module, total height: 1650 mm, see sketch of oven dimensions G/H/I) Width/Height/Depth			529/450/ 383 V 820	529/290/ 463 2020	689/13 533 2420	
5 L. L.		Electrical load (maximally equipped) at 230 V, 50/60 Hz approx. W					
Packing data		Net weight approx. kg			83 104	110 135	
	Gross weight (packed in carton) approx. kg					830/10	
	Packed dimensions (Width, Height, Depth) approx. mm					800 41/5	
	Net weight pump module without/with Gross weight pump module without/wi	· ·	approx. k	-	25/41 30/46 46/62 51/67		
			approx. kg		660/870/590		
Packed dimensions pump module (Width, Height, Depth) approx. mm					000/8/0/590		
Order No. Vacuum Ovens						V010	
ptions			29	49		101	
Premium module: comprises the inert gas inlet (only size 49 and 101), extra connectors for thermoshelves, 1 (size 29), 2 sizes 49/101)					T5		
- 20 mA current	- 20 mA current loop interface (requires option T5)  Temperature actual value (0 to 210 °C = 4 - 20 mA)				V3		
Temperature of a Pt100 sensor positioned flexibly in chamber for external temperature monitoring - price per sensor				V6	V6		
Potential-free contact (24 V/2 A) with socket, for combination error message (e.g. supply failure, sensor fault, fuse)					H6		
MobileALERT, notification by SMS in case of any error or alarm of the device (requires option H6)				C3			
nobileALERI for 2 ption T5)	? alarm notifications; notification by SMS. ten	nperature and vacuum alarm (only in connection with	-		C4		
Vorks calibration ustomer specifica	certificate for one (freely selectable) temperation	ture and pressure value (per thermoshelf) according to		D00116			
ccessories				29	49	10	
Fhermoshelf - aluminium eloxadised material WSt. 3.3547 (ASTM B209) with integrated large-area heating including local temperature sensing Pt100, 4-wire-circuit); individual overtemp. protection for each shelf MLOP (Multi-Level-Overtemperature-Control) and calibration certificate					B00743	B007	
Thermoshelf - stainless steel material 1.4404 (ASTM 316 L) for especially corrosive material with integrated large-area heating including local temperature sensing (Pt100, 4-wire-circuit); individual overtemp. protection for each shelf MLOP (Multi-Level-Overtemperature-Control) and calibration certificate					B00734	B007	
Subframe, tubular steel, black enamelled (for stacking unit consisting of vacuum oven and pump module, total height: 1650 mm, see "further data" and sketch of oven dimensions)					E02031	E020	
and sketch of oven dimensions) Works calibration certificate for 3 temperatures: +50, +100, +160 °C at 20 mbar pressure. Price per thermoshelf					D00115		
Guarantee extension by 1 year					GA2Q5		
Noise-insulated vacuum pump module without pump (exterior dimensions and -material No. s. vacuum oven) with antivibration metal plate at the bottom to accommodate the vacuum pump, incl. full-sight glass door. Socket, signal cable and connecting hose to the vacuum oven					PM49	PM1	
bottom to accommodate the vacuum pump, incl. full-sight glass door. Socket, sighal cable and connecting hose to the vacuum oven  Noise-insulated vacuum pump module, as above, however with built-in pump, 230 V, 50/60 Hz					PMP49	PMP1	
Signal cable (3 m) for control of rotation speed and optimising pump performance by demand-controlled activation of purge of Memmert pump (not required with pump module)					PMP29   PMP49   PM B39410		
acuum connectir ump module)	ng hose (3 m) from oven to Memmert pump	incl. optimised connection accessories (partially stainless stee	el), (not required with		B04026		
Vacuum pump with chemically resistant 4x diaphragm, pump capacity at atm. pressures: approx. 50 NI./min = 3,0 m <sup>3</sup> /h and autom. purge control, 230 V, 50 Hz. Max. guarantee period 2 years (requires acessories B39410 and B04026)					E07509		
USB-Ethernet adapter					E06192		
Ethernet connection cable 5 m for computer interface  USB User-ID stick (with User-ID licence): Oven-linked authorisation licence (User-ID-programme) on Memory-stick, prevents undesired manipulation					E06189		
y unauthorised t	hird parties. When reordering please specify	serial number		1011	B33170		
FDA conforming software AtmoCONTROL (FDA edition). Meets the requirements for the use of electronically stored data sets and electronic signatures as laid down in Regulation 21 CFR Part 11 of the US Food and Drug Administration (FDA). Base licence for the control of one unit. Respective IQ/OQ documents available in German and English language (without surcharge)					FDAQ1		
Integration of additional units (up to max. 31 units) into an already existent FDA-software licence					FDAQ2		
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer					D00124		
IQ/OQ document with device-specific works test data for one free-selectable temperature and vacuum value, incl. temperature distribution survey at Memmert for 5 measuring points per thermoshelf to DIN 12880:2007-05. PQ check list as support for validation by customer valid for one thermoshelf; dito further thermoshelves, price on demand. 360 € for further temperature and vacuum values					D00117		
On-site IQ/OQ for a freely selectable temperature and pressure value of a thermoshelf, including temperature distribution survey for 5 measuring points to DIN 12880: 2007-05 (excluding travel costs, not subject to discount, GER, AT, FR only)					DLQ106		
Extension of DLQ106 by an additional freely selectable temperature and pressure value (each thermoshelf) (not subject to discount)					DLQ106A		
Extension of DLQ106 by a further thermoshelf for a freely selectable temperature and pressure value (not subject to discount)					DLQ106T		
Individual on-site Performance Qualification (PQ)  Maintenance VO - carrying out and documentation according to Memmert maintenance plan (excluding travel costs, not subject to discount, GER,					DLQ200 S00320		
AT, FR only) Maintenance cont	tract VO - carrying out and documentation ac	cording to Memmert maintenance plan, minimum duration	3 years (excluding trav	el	S00320J		
costs, not subject to discount, GER, AT, FR only)  Calibration of one freely selectable temperature value (excluding travel costs, not subject to discount, GER, AT, FR only)					S00205		
Calibration of an additional temperature value (not subject to discount)					S00215		
			FR only)		500212		